



Exporting Setting and/or history to USB

1. When the system is ready, enter the QUICK OPTIONS screen by pressing the UP and DOWN keys simultaneously.

You will see the quick options screen.

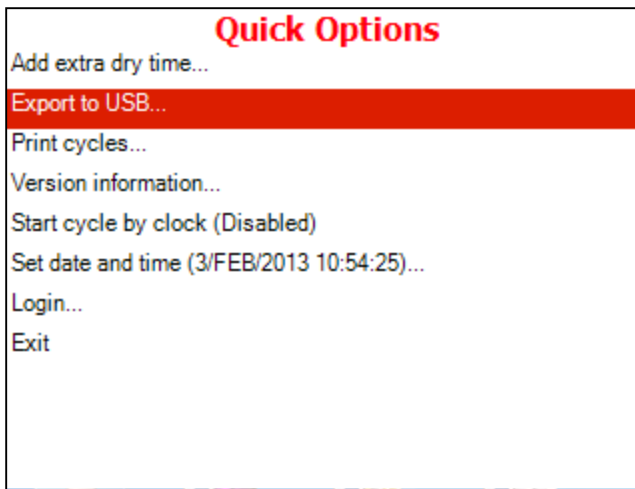

2. Move the cursor to Export to USB
3. Press the START/STOP key.
4. The following screen will be displayed:

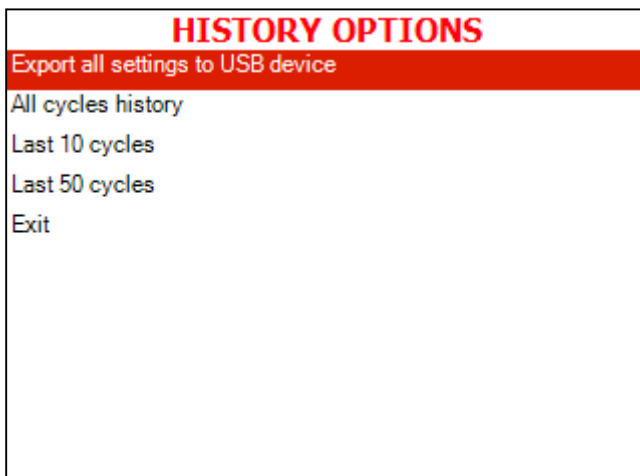

These options enable to export of both the history and settings to the USB device.


Please note that the settings and the history exports are two separate exports so if you need both groups you will need to do two exports. The settings export will generate two folders, "FactoryDefault" and "Current", while the history export will generate a "History" folder.

5. Insert the USB device into the USB Socket
6. With the up/down arrow key, move the cursor to the required item (either "Export all settings" or one of the history exports) and press the **START/STOP** key.
7. The following screen will be displayed:




8. If another export is required, repeat the above steps and select the appropriate export
9. When finished, remove the USB device from the USB Socket.
10. In order to exit this subdirectory, press the START/STOP key.


In order to email these files to us. Go into your USB key, select all the folders that have been created by the download from the autoclave.

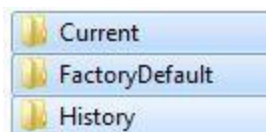

Once they have been selected you should be able to right-click on them and send them to a compressed/ZIP file.

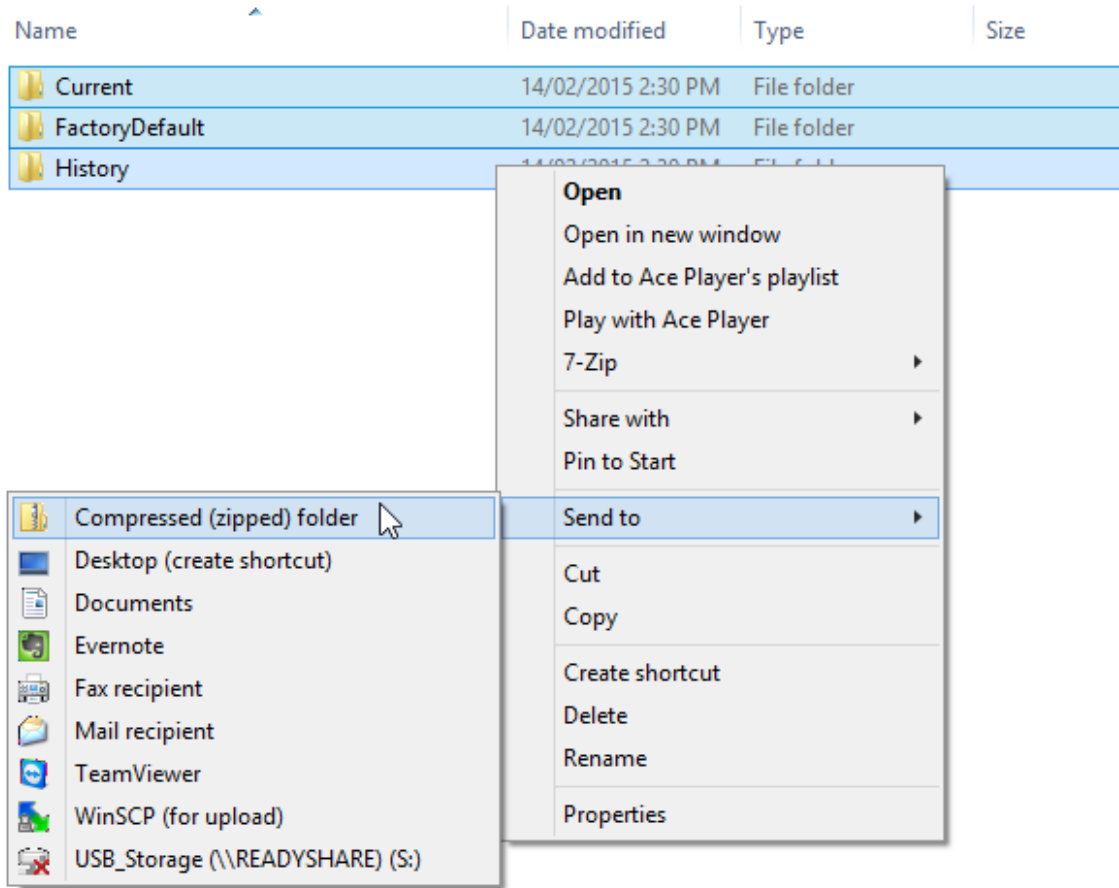

Once you have done this you will end-up with a single compressed .ZIP file that you can email to us.

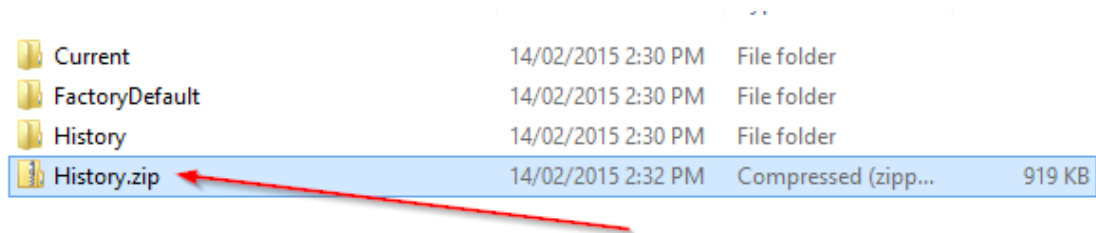

The actual name of the .zip file may vary depending on the order of your folders. Please email this file to us at service@labgearaustralia.com.au or the address provided.