


Import Application From USB

Unzip the supplied application folder compressed file

Use a clean USB key and copy the unzipped application folder to the root directory of the USB key

Insert the USB key into the autoclave

Log in to Technician mode (code "0321") and go to Version Handling

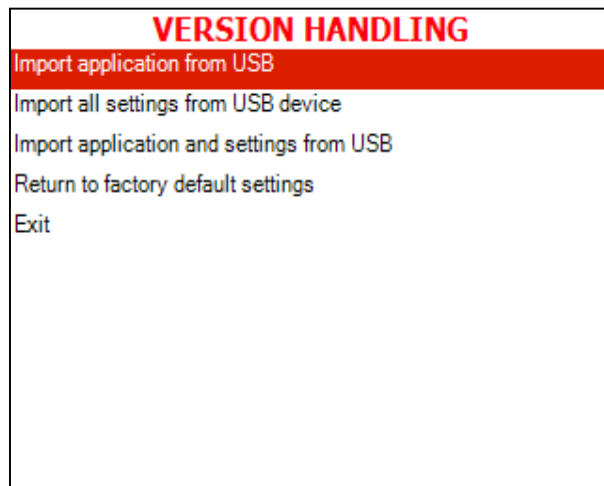

Import application from USB

In this menu you can replace the autoclave software with the application software from the USB device to the autoclave.

1. Browse to the menu item:

Version handling\Import application from USB

The system will prompt you to confirm import

2. Move the cursor to Confirm and press Start/Stop.

The application will be imported from the USB device.