

SNIJDERS EVOSAFE FREEZER USER SETTINGS ADJUSTMENT

To View Temperature Set Point	Press & Hold SETP
To Mute Alarm	Press Mute/Battery
How To Enter Access Code	Press Code Key
	Enter Code 1234 (+, SETP, MUTE, -)
	Wait for display to stop blinking. Code light will remain lit & you are now in the user level access area
To Clear Error	Enter Access Code (1234)
	Press SETP for 2 seconds
	<i>(note if error still valid it will not clear)</i>
To Change Setpoint	Enter Access Code (1234)
	Press and hold SETP
	Adjust setpoint via +/- keys
To Adjust Cold Alarm Setpoint	Enter Access Code (1234)
	Enter Parameter Mode by pressing PRG Buton for 2 seconds
	Parameter P30 will show on display
	Press and hold SETP
	Adjust setpoint via +/- keys
To Adjust Warm Alarm Setpoint	Enter Access Code (1234)
	Enter Paramater Mode by pressing PRG Buton for 2 seconds
	Parameter P30 will show on display
	Go to P31 via +/- keys
	Press and hold SETP
	Adjust setpoint via +/- keys
To Access Parameters for Adjustment	Enter Access Code (1234)
	Press Program Button for 2 seconds
	Use +/- key to scroll through parameter numbers
	Hold SETP Button to view parameter
	Hold SETP Button then +/- to adjust parameter
USER ADJUSTABLE PARAMETERS	
P30 - Cold Alarm Set Point	P50 - CO2/LN2 Set Point
P31 - Warm Alarm Set Point	P70 - Alarm Relay Delay
Alarm Codes	
Hi 1 - Warm Temperature Alarm	E5 - Battery Voltage Low
Hi 2 - Condensor Temperature too warm	E6 - Door Alarm
Lo 1 - Cold Temperature Alarm	E10 -Mains voltage out of range
USB Download	Insert USB Key in port behind condensor panel
	Enter Access Code (1234)
	Press the + key
	Wait for L.Usb to appear on the screen
	Press SETP
	Screen should now count up to 100% for download progress
	Once download complete remove USB Key